

Pulseless Dosing Disposable Pump compatible with Cole Parmer - Masterflex L/S drives.



www.naopump.com

User manual

1. Position the holder

2. Fix the 4x screws

3. Insert the Rotor


(Do not remove the Materflex motor shaft rubber)




4. Set the Masterflex Driver parameters

| | | |
|----------------|----------------------|----------------------------------|
| Tubing size | 36 | |
| Calibration | 333.333 | Max. 100 RPM → 0.5 to 500 ml/min |
| Pump direction | CW only (Clock Wise) | |

5. Insert the Disposable and lock the red door

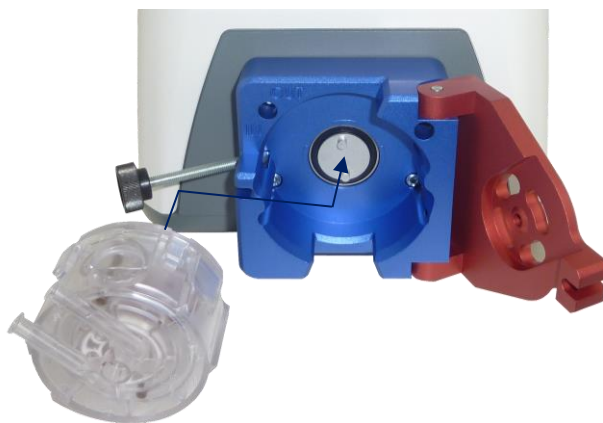

6. Connect the inlet and outlet tubing (Luer Lock Male)

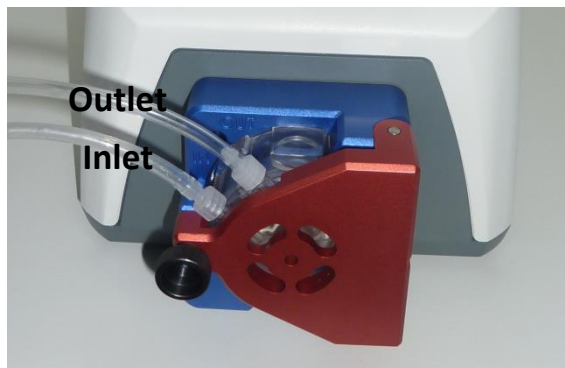

The pump is ready